


Easy Opening of Push-Tear-Off Caps


1/ Lift the plastic lid with your thumb at the triangle until the lid is in a vertical position

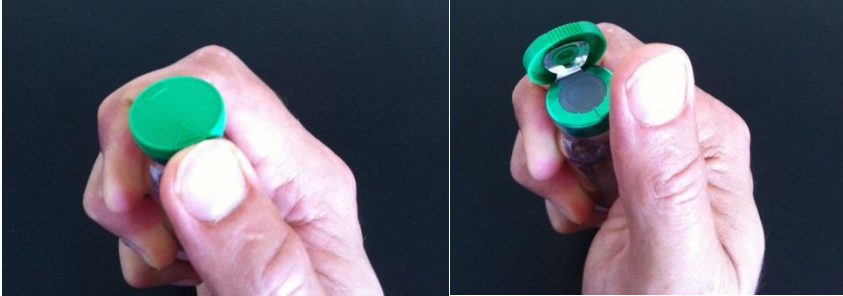

2/ Tear straight downwards until the right aluminum-attach opens



3/ Tear to the left (in direction of the arrow) to remove the aluminum cap completely

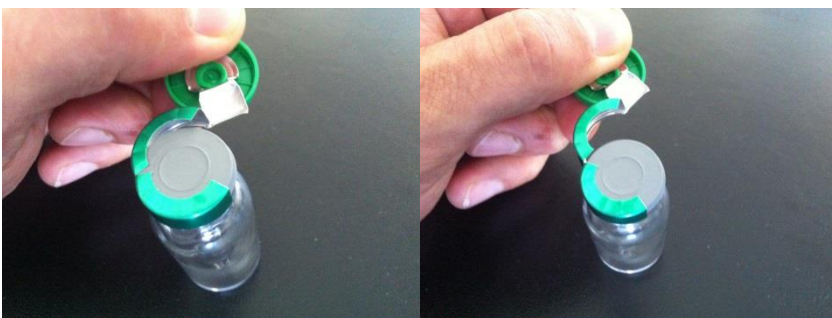

EMA Pharmaceuticals S.A.S.

460 Route de Monçay - 45740 LAILLY EN VAL - FRANCE

Tél : + 33 2 38 46 96 60 Fax : + 33 2 38 44 70 88

contact@emapharma.com

www.emapharma.com